

# Prizmatix

## UHP-Microscope-LED-435 Ultra High Power LED Light Source for Fluorescence Microscopy

Ver. 02

### Introduction

The Ultra High Power LED (29 Watt) provides excellent replacement of conventional arc lamp sources in fluorescence microscopy applications like Cyan GFP (CFP) stain. Many power demanding applications such as chemical reaction activation, machine vision, and numerous others may benefit of this new LED light source. The UHP-Mic-LED-435 provides **>1Watt of collimated** LED power on its output. The LED driver supports CW or pulsed operation.



### Key Features

- Compatible with Prizmatix modular Microscope-LED Light Source products family
- Single chip Ultra High Power LED
- TTL external triggering
- Excellent for fluorescence excitation
- LED spectrum can be narrowed by optional band pass filter
- Stable precisely adjustable power
- Long life (no lamp or laser tube replacement required)
- Speckle free
- Rapid warm up time

### Applications

- Fluorescence microscopy with fluorophores CFP (cyan GFP), ECFP, AmCyan, SYTOX Blue, BOBO-1, BO-PRO-1, DEAC, Quantum Dots
- Whole body imaging of small animals in-vivo
- Photo activation (PA)
- Fluorescence recovery after photo bleach (FRAP)
- Fluorescence in situ hybridization (FISH)
- Bio analysis
- Machine Vision

#### Main Office

Phone: +972-72-2500097  
Fax: +972-72-2500096  
sales@prizmatix.com

#### European Sales Office

Phone: +44 (0)77-9172-9592  
Fax: +44 (0)20-7681-2977  
sales.europe@prizmatix.com

#### North America Sales Office

Phone: +1-(248)-436-8085  
Fax: +1-(248)-281-5236  
sales.usa@prizmatix.com

P.O.B. 4234 Modiin-Ilite 71919, Israel

# Prizmatix

## Optical Specifications

Wavelength	nm	435
Wavelength range	nm	±5
Spectrum half width	nm	16.5
Collimated optical power output	mW	>1000

## Benchtop LED Current Controller Specifications

### Features

- Constant current or chopping modes
- Precise LED current setting by 10 turn dial
- TTL external trigger input
- Compact and robust enclosure

Digital modulation input		TTL
Connector for TTL input		BNC
Digital modulation frequency	Hz	DC-30000
Rise / Fall time (10% - 90%)	µs	2
ON/OFF extinction		>10 <sup>6</sup> : 1
Input Voltage	V	12
Max Input current	A	5.5
Power Adaptor Input		100-240 VAC, 50-60Hz, 1.5A

## General Specifications

Operation temperature range	°C	10-35
Storage temperature range	°C	-10-55
Operating relative humidity (Non condensing)	%	<90
Head dimensions		See drawing below
Head weight	g	500
Controller dimensions (LxWxH)		220mm x 165mm x 51.5mm
Controller weight	g	700
Power adaptor dimensions (LxWxH)		146mm x 76mm x 44mm
Power adaptor weight	g	600
Fan noise	dBA	38

### Main Office

Phone: +972-72-2500097  
Fax: +972-72-2500096  
sales@prizmatix.com

### European Sales Office

Phone: +44 (0)77-9172-9592  
Fax: +44 (0)20-7681-2977  
sales.europe@prizmatix.com


### North America Sales Office

Phone: +1-(248)-436-8085  
Fax: +1-(248)-281-5236  
sales.usa@prizmatix.com

P.O.B. 4234 Modiin-Ilite 71919, Israel


# Prizmatix

## Performance

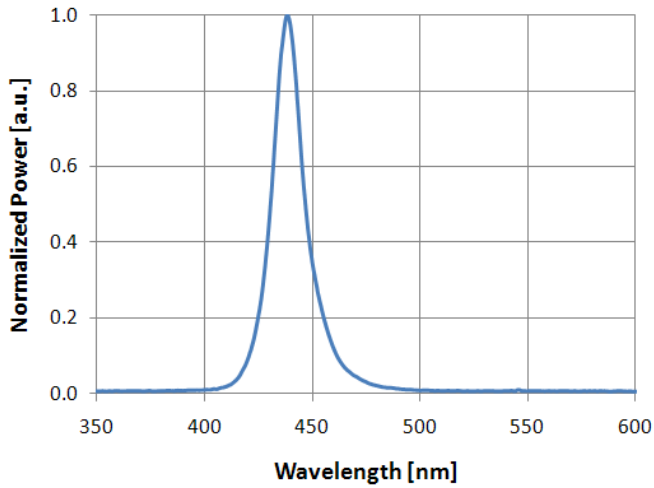

- UHP-Microscope-LED-435 spectrum

## Mechanical Drawings

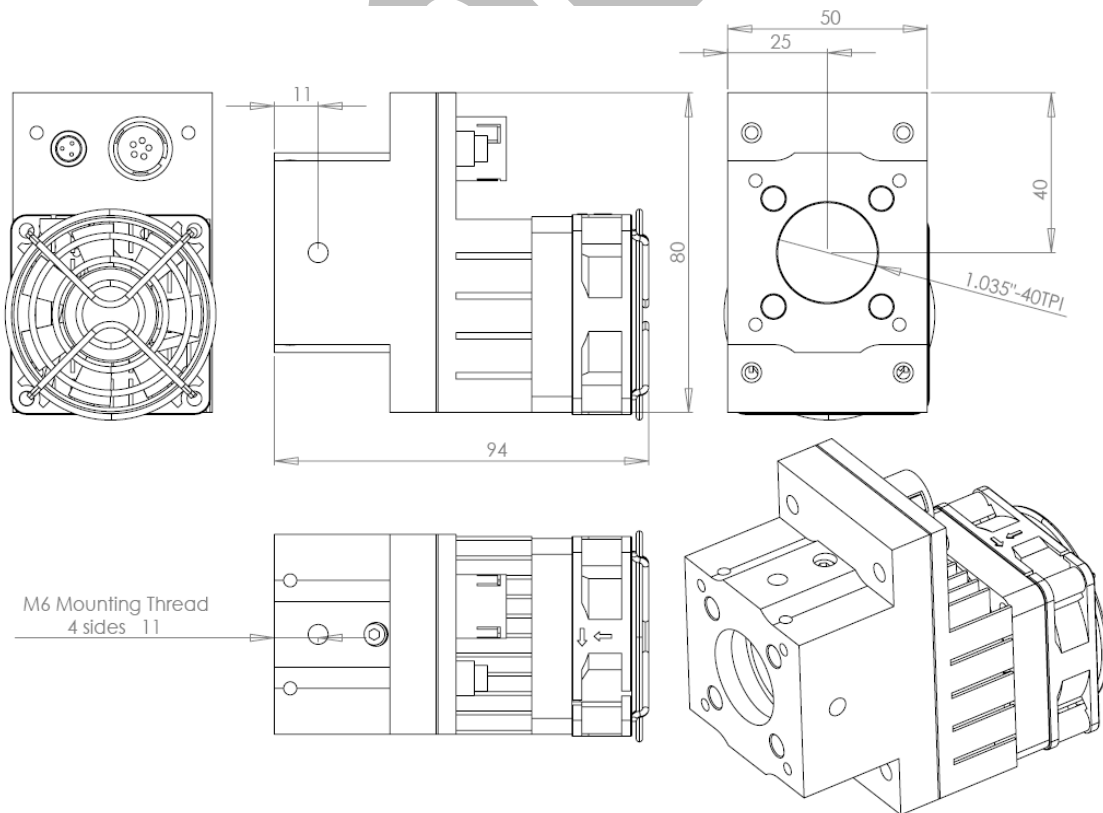

### Main Office

Phone: +972-72-2500097  
Fax: +972-72-2500096  
sales@prizmatix.com

### European Sales Office

Phone: +44 (0)77-9172-9592  
Fax: +44 (0)20-7681-2977  
sales.europe@prizmatix.com

### North America Sales Office

Phone: +1-(248)-436-8085  
Fax: +1-(248)-281-5236  
sales.usa@prizmatix.com

P.O.B. 4234 Modiin-Ilite 71919, Israel

# Prizmatix

## Optional Accessories

### Filter Wheel:

The UHP-Mic-LED can be equipped with a 6 positions filter wheel at the beam output. This accessory is especially useful for UHP-Mic-LED-White light source. Please see video clip <http://www.youtube.com/watch?v=iv7dlwLHaUE> for more details.

### Fiber Coupler Adaptor:

The UHP-Mic-LED can be easily changed from direct microscope coupling to fiber coupled LED configuration by means of Fiber Coupler Adaptor (SMA, CF or ST connector). This adaptor can be easily assembled by a user on the Mic-LED output. Please see video clip <http://www.youtube.com/watch?v=iv7dlwLHaUE> for more details.

### Liquid Light Guide Adaptor:

The Microscope-LED can be easily changed from direct microscope coupling to Liquid Light Guide coupled LED configuration by means of LLGA Adaptor. This adaptor can be easily assembled by a user on the Mic-LED output. Please see video clip <http://www.youtube.com/watch?v=iv7dlwLHaUE> for more details.


### Collimator:

The output from optical fiber is divergent according to fiber NA. In order to reduce the divergence angle a collimator module can be used. Prizmatix collimator was especially designed to fit thick core high NA Polymer Optical Fibers. See more details at <http://prizmatix.com/docs/collimator.cfm>

### Fiber Bundles:

To combine outputs of multiple LEDs a Y-shaped fiber bundle with two or more input branches can be used. Prizmatix can help to configure and build custom fiber bundles for specific applications. See more details at:

[http://prizmatix.com/docs/Custom\\_Fiber\\_Optic\\_Assemblies.cfm](http://prizmatix.com/docs/Custom_Fiber_Optic_Assemblies.cfm)

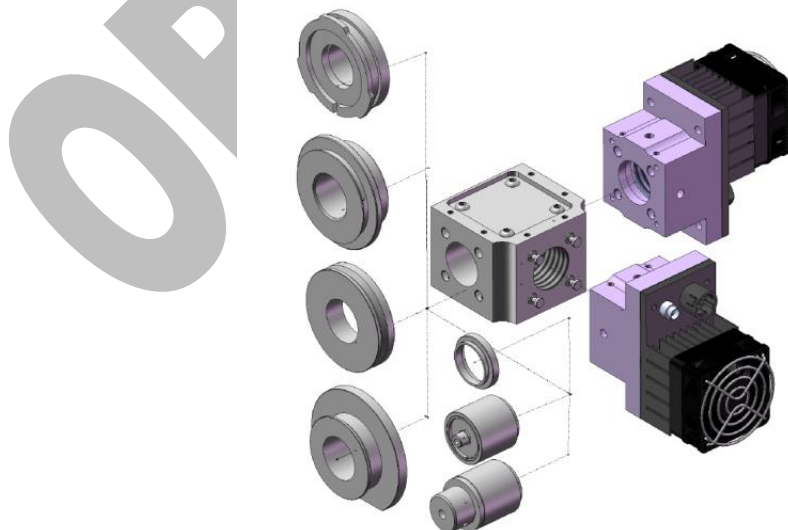

---

### Main Office

Phone: +972-72-2500097  
Fax: +972-72-2500096  
[sales@prizmatix.com](mailto:sales@prizmatix.com)

### European Sales Office

Phone: +44 (0)77-9172-9592  
Fax: +44 (0)20-7681-2977  
[sales.europe@prizmatix.com](mailto:sales.europe@prizmatix.com)

### North America Sales Office

Phone: +1-(248)-436-8085  
Fax: +1-(248)-281-5236  
[sales.usa@prizmatix.com](mailto:sales.usa@prizmatix.com)

---

P.O.B. 4234 Modiin-Ilite 71919, Israel

---